

PIG Biohazard Spill Kit

Instructions

Step 1 – Wear appropriate PPE

Put on the personal protective equipment supplied in the spill kit before cleaning up any fluids.

Step 2 – Clean up the spill

For small spills, slowly pour the PIG Absorb-&-Lock Biofluid Absorbent onto the spill area, working from the outside inward until it's completely absorbed and solidified in 2–3 minutes. Remove with scraper and pan (included).

For larger spills, use the PIG Absorb-&-Lock Biofluid Absorbent and PIG HazMat Pads to absorb the spilled liquids.

Step 3 – Sanitize the surface

Use the premoistened surface sanitizing wiper to remove any residue from the surface.

Step 4 – Dispose of cleanup materials

Place all used articles in the red biohazard disposal bag and tie the bag closed. (See disposal notice below.)

Step 5 – Sanitize hands

Use the premoistened hand sanitizing wiper following removal of PPE.

CAUTION: Assume all blood and body fluids are contaminated and potentially hazardous to your health.

NOTE: Store product in a cool, dry and well-ventilated area.

DISPOSAL NOTICE: Used absorbents can be considered hazardous and must be disposed of in compliance with local, state and federal regulations. For information on disposal options, call New Pig Technical Services at 1-800-HOT-HOGS (468-4647).



New Pig

One Pork Avenue
Tipton, PA 16684-0304

newpig.com

North America: **1-800-468-4647**

Europe: **+31 (0)76 596 92 50**

PIG, PIG logo are registered trademarks in USA and other countries. See tm.newpig.com

UK: **0800 919 900**

Outside North America: **+1-814-684-0101**

165916 / 331009