

# PIG BioSafety Spill Cleanup Kit

## Instructions

### Step 1 – Wear appropriate PPE

Put on the personal protective equipment supplied in the spill kit before cleaning up any fluids.

### Step 2 – Clean up the spill

**For small spills**, slowly pour the Red Z Absorbent onto the spill area, working from the outside inward until it's completely covered. The liquid will be completely absorbed and solidified in 2–3 minutes. Remove with scraper and pan (included).

**For larger spills**, use the PIG Haz-Mat Pads to absorb the spilled liquids. Make a final pass with dry wiping cloth (included).

### Step 3 – Sanitize the surface

Use the pre-moistened surface sanitizing wiper to remove any residue from the surface.

### Step 4 – Dispose of cleanup materials

Place all used articles in the red biohazard disposal bag and tie the bag closed. (See disposal notice below.)

**CAUTION:** Assume all blood and body fluids are contaminated and potentially hazardous to your health.

**NOTE:** Store product in a cool, dry and well-ventilated area.

**DISPOSAL NOTICE:** Used absorbents can be considered hazardous and must be disposed of in compliance with local, state and federal regulations. For information on disposal options, please call New Pig Technical Services at 1-800-HOT-HOGS (468-4647).



Specialized kit includes PIG Absorbents, blood solidifier and other tools to help you quickly and safely clean up biofluids.



## New Pig

One Pork Avenue  
Tipton, PA 16684-0304

**Still have questions? In North America contact us:**

By Phone: **1-800-HOT-HOGS** (468-4647)

Online: **newpig.com**

Email: **hothogs@newpig.com**